

Method Procedure



Document Number: SASoM/METHOD/067.v5

Title: Preparation of Saturated Bromophenol Blue Solution

Version: v5

Author: Peter Mullen

Effective from:	11/11/2022	
Valid to:	11/11/2024	

SOP History		
Number	Date	Reason for Change
v1	15/07/2014	Original
V2	15/07/2016	Biennial Review
V3	15/07/2018	Biennial Review
V4	11/11/2020	Biennial Review
V5	11/11/2022	Biennial Review

1.0 Purpose -

This SOP describes the current procedure for preparing saturated bromophenol blue solution in Laboratory 248 at the St Andrews School of Medicine (SASoM).

2.0 Scope -

This SOP applies to the staff in the SASoM involved in preparing saturated bromophenol blue in Lab 248.

3.0 Responsibilities -

All staff involved in cell culture are responsible for ensuring that the methods are followed in accordance with this SOP.

All staff must have read and signed the relevant risk assessment documents before performing this procedure.

St Andrews School of Medicine (SASoM) Systems Pathology Group

Method Procedure



4.0 Procedure -

- 1. Add a heaped spatula full of solid bromophenol blue powder to distilled water in a universal container.
- 2. Mix
- 3. Continue to add more bromophenol blue.
- 4. Stop adding when no more will dissolve.
- 5. Centrifuge to pellet the undissolved powder.
- 6. Carefully transfer the saturated supernatant to a fresh tube.

5.0 Personal protection -


A Howie laboratory coat and lab gloves must be worn at all times.

6.0 Spillages -

Always clean up any spills immediately after use, only you know what you have spilt and are aware of its hazard. Spillages should be mopped up with paper towel, disinfected with 70% ethanol and finally washed with warm water.

7.0 Training -

8.0 Related documents - RA28951 SDS-PAGE Western Blotting.

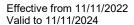



St Andrews School of Medicine (SASoM) Systems Pathology Group

SCHOOL OF MEDICINE

Method Procedure

9.0 Approval and sign off -

Author:

Name: Peter Mullen

Position: Research Fellow

Signature: Notes Muller

Date: 28/11/2022

Management Approval:

Name: David Harrison

Position: Professor

Signature: Date: 25/01/2023

QA release by:

Name: Peter Mullen

Position: QA Manager

Signature: Voter Wuller Date: 28/11/2022

St Andrews School of Medicine (SASoM) Systems Pathology Group



Method Procedure

STANDARD OPERATING PROCEDURE

Please sign below to indicate you have read this S.O.P and understand the procedures involved.

NAME	POSITION HELD	SIGNATURE	DATE